




Analyses Report

Determination of Nitrogen on Dog Crunches

*R&D Department
NC Technologies S.r.l.*




Typical calibration curve with certified standard

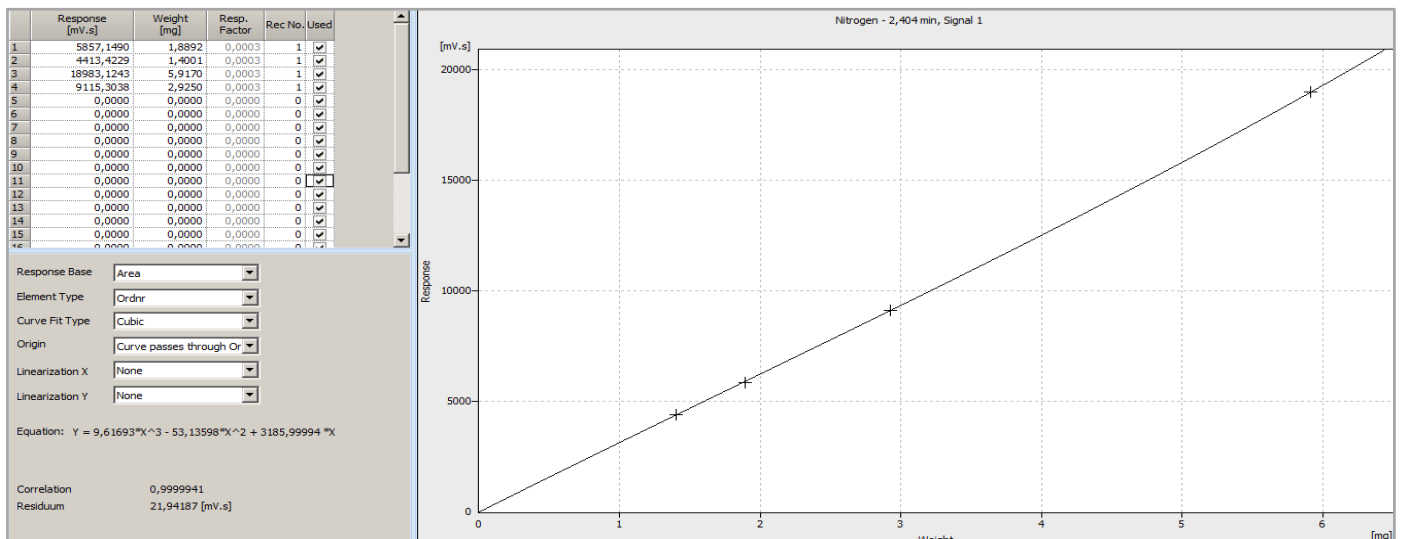

Calibration curve for N; range 1mg to 6mg
Standard: Acetanilide


EDTA vs Dog Crunches: verification of instrument and method

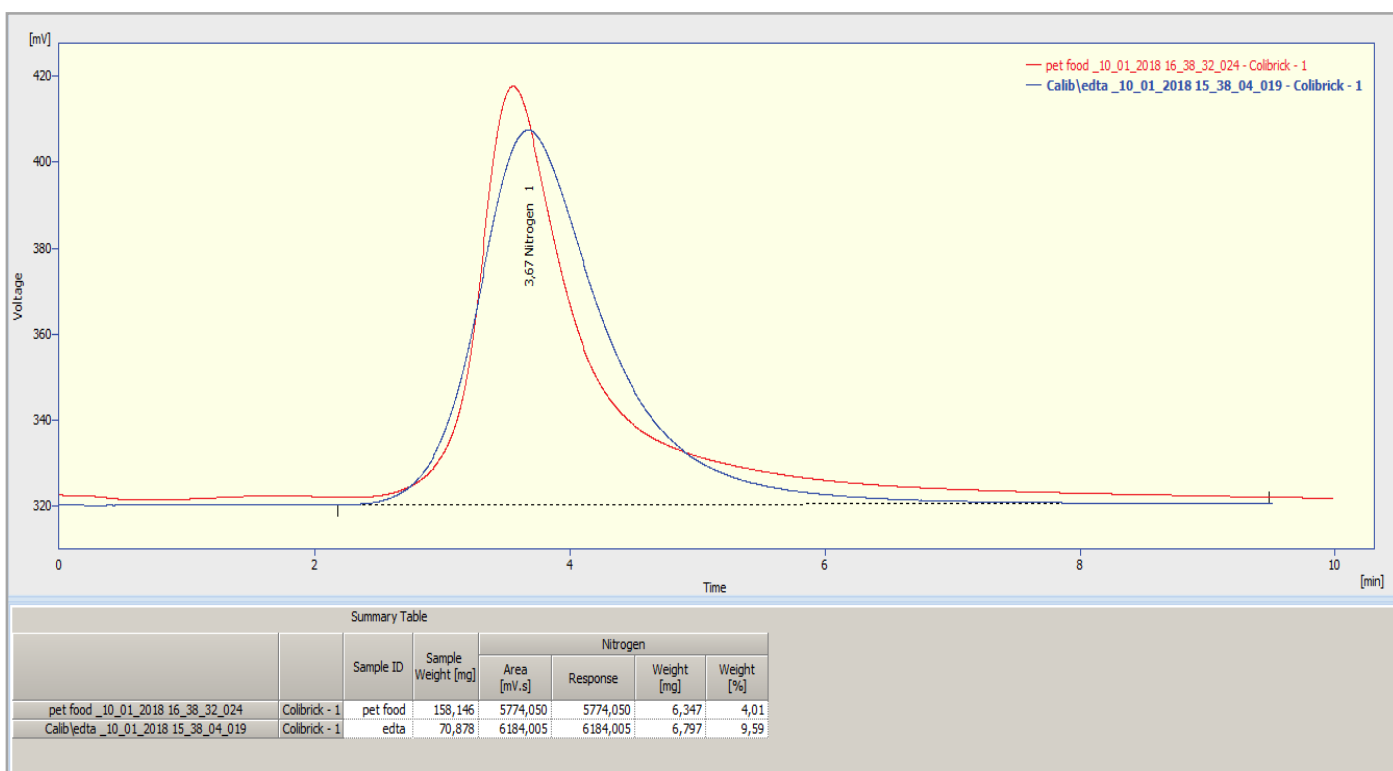

Possibility of high sample dosage with the same analysis performances



Dog Crunches nitrogen determination: different weights

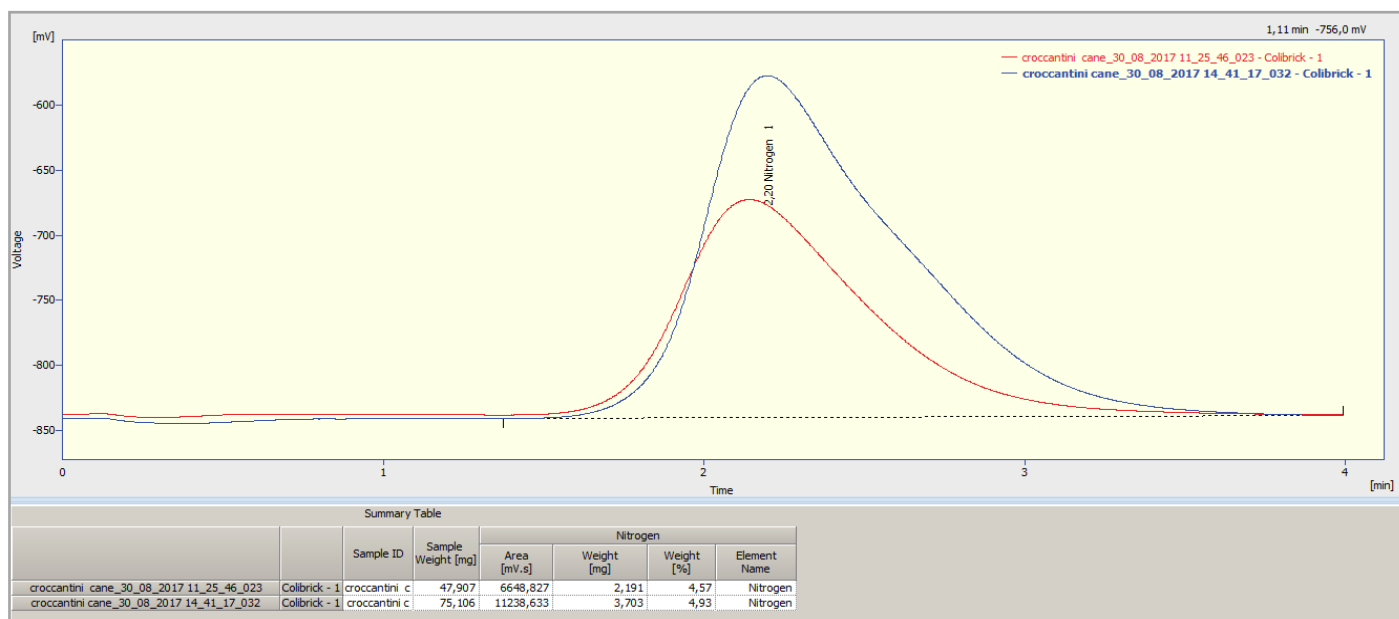

Dog crunches	Sample weight [mg]	Nitrogen Weight [mg]	Nitrogen Weight [%]
	47.907	2.191	4.57
	75.106	3.703	4.93